



# **CITRAD™**

## **Acid Cleaner & Detergent**

---

### **- BRIEF DESCRIPTION -**

*Citrad concentrated acidic detergent is ideal for the removal of salts, scale, and metal-oxide. Also used for metal brightening and pharmaceutical-process equipment. Citrad contains a safe, organic acid and is phosphate-free, so it will not interfere with sensitive equipment or testing procedures. Citrad is also completely free-rinsing and biodegradable.*

### **- USE PROTOCOL -**

Citrad concentrate must be diluted with distilled, deionized or clean potable water and can be used as a soak, scrub or ultrasonic cleaner.

Citrad is used to clean labware, stills, cleanrooms, optical parts, pharmaceutical equipment, and manufacturing equipment. Citrad is also effective in removing metal oxides, scale, soil, grease, chemicals and solvents and other compounds from surfaces and equipment.

### **- DIRECTIONS FOR USE -**

Add 1.25 ounces to 2.5 ounces per gallon of water (1% to 2% v/v). For more difficult applications increase the concentration of Citrad solution and/or increase the temperature.

Citrad will foam if dispensed by sprayer.

### **- DISPOSAL -**

All ingredients in Citrad are biodegradable and are considered safe when recommended use dilutions are disposed of by drain.

### **- PACKAGING -**

**1 Gallon / 4 per case**

### **- CATALOG NUMBER -**

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	1 Gal.	4401

---

**For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

---

®Decon Labs, Inc.

Revised 8/18/2011