



Contrex® AL is a concentrated alkaline liquid detergent. It removes virtually any contaminant including grease, oils, blood, tissue, particulates, deposits, chemicals and solvents. Contrex® AL works by soaking, scrubbing, or in ultrasonic cleaning baths. Contrex AL is totally rinsable eliminating detergent build-up that interferes with analysis, tissue culture, and other critical procedures. Its activity can be accelerated by heating, agitation, and ultrasonics, or by increasing concentrations. Contrex® AL is phosphate-free, biodegradable and readily disposable.

Packaging		
Size	Case QTY	Decon #
1 Quart	12	5332
1 Gallon	4	5301

APPLICATIONS: Contrex® AL is not to be used with machines, product will foam. After cleaning, rinse thoroughly. Safe on most surfaces including glass, plastic, and most metals. Cleaning applications include:

- Labware
- Instruments
- Equipment
- Production Mixers/fillers/storage containers
- Optics

STORAGE: Keep contents in a clean and dry environment.

WORKING SOLUTIONS

Make a fresh 1% Contrex® AL working solution (1.25oz. per gallon) in deionized or distilled water for best results. Soak, scrub or sonicate as needed.

RINSABILITY

Contrex® AL is totally rinsable - however, numerous factors affect rinsability. These factors include water quality, temperature, volume of water, nature of surface, shape/size of the vessel.



For details on Decon's Full Line of Detergents visit

www.deconlabs.com

Rev. 09/2022