



Technical Data Sheet

- BRIEF DESCRIPTION -

Contrex AL is a concentrated alkaline liquid detergent. It removes virtually any contaminant including grease, oils, blood, tissue, particulates, deposits, chemicals and solvents. Contrex AL works by soaking, scrubbing, or in ultrasonic cleaning baths. Contrex AL is totally rinsable eliminating detergent build-up that interferes with analysis, tissue culture, and other critical procedures. Its activity can be accelerated by heating, agitation, and ultrasonics, or by increasing concentrations. Contrex AL is phosphate-free, biodegradable and readily disposable.

- APPLICATIONS -

Not for use with machines, product will foam. After cleaning rinse thoroughly. Safe on most surfaces including glass, plastic, and most metals. Cleaning applications include:

- Labware
- Instruments
- Process Equipment
- Production mixers, fillers, storage containers
- Optics

- WORKING SOLUTIONS -

Make up a fresh 1 % Contrex AL working solution (1.25 oz per gallon) in deionized or distilled water for best results. Soak, scrub or sonicate as needed.

- RINSABILITY -

Contrex AL is totally rinsable – however, numerous factors affect rinsability. These factors include temperature, and volume of water, nature of surface, shape and size of the vessel.

-STORAGE -

Product must be stored in a clean, dry environment.

- PRECAUTIONARY STATEMENTS -

KEEP OUT OF REACH OF CHILDREN CAUTION

Contents may cause eye irritation. Avoid contact with eyes.

-FIRST AID-

If in eyes: flush eyes with copious amounts of water for 15 minutes. Seek medical attention if eye irritation persists.

- PACKAGING -

1 Quart / 12 per case
1 Gallon / 4 per case

- CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
1 Quart	5332
1 Gallon	5301

For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or

See our website at www.deconlabs.com

©Decon Labs, Inc.

Revised 4/02/2012