



Technical Data Sheet

- BRIEF DESCRIPTION -

Contrex AP is a powdered labware detergent. It removes virtually any contaminant including grease, oils, blood, tissue, particulates, deposits, chemicals and solvents. Contrex AP works by soaking, scrubbing, or in ultrasonic cleaning baths. Contrex AP is totally rinsable eliminating detergent build-up that interferes with analysis, tissue culture, and other critical procedures. Its activity can be accelerated by heating, agitation, and ultrasonics, or by increasing concentrations. Contrex AP is biodegradable and readily disposable.

- APPLICATIONS -

Not for use with machines, product will foam. After cleaning, rinse thoroughly. Safe on most surfaces including glass, plastic and most metals. Cleaning applications include:

- Labware
- Instruments
- Process Equipment
- Production mixers, fillers, storage containers
- Optics

- WORKING SOLUTIONS -

Make up a fresh 1 % Contrex AP working solution (1.25 oz per gallon) in deionized or distilled water for best results. Soak, scrub or sonicate as needed.

- RINSABILITY -

Contrex AP is totally rinsable – however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape and size of the vessel.

- STORAGE -

Keep contents in a clean and dry environment.

- PACKAGING -

4 lb. /9 per case
25 lb. /each
50 lb. /each

- CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
4 lb.	5204
25 lb.	5225
50 lb.	5250

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

®Decon Labs, Inc.

Revised 4/02/2012