



Technical Data Sheet

- BRIEF DESCRIPTION -

Contrex CA concentrated acidic detergent is formulated for use in descaling, mineral removal, oxide removal or in applications where an acidic detergent is indicated. Contrex CA contains a safe, organic acid and is phosphate-free, so it will not interfere with sensitive equipment or testing procedures. Contrex CA is also completely free-rinsing and biodegradable.

- APPLICATIONS -

Contrex CA concentrate can be diluted with distilled, deionized or clean potable water and can be used as a soak, scrub or ultrasonic cleaner.

Contrex CA is not for use in machines-product will foam. In order to accelerate cleaning increase the solution strength or temperature. Contrex CA is safe on most surfaces including glass, plastic and most metals.

Cleaning applications include:

- Labware
- Trace metal, scale and metal-oxide removal
- Metal brightening
- Process equipment
- Optical
- Scale removal

- WORKING SOLUTIONS -

Make up a fresh 1 to 2 % Contrex CA working solution (1.25 to 2.5 oz. per gallon) in deionized or distilled water for best results. Clean by scrubbing, soaking or by ultrasonic cleaning. For critical applications use deionized or distilled water. After cleaning, rinse thoroughly.

- RINSABILITY -

Contrex CA is totally rinsable – however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape and size of the vessel.

- DISPOSAL -

All ingredients in Contrex CA are biodegradable and are considered safe when recommended use-dilutions are disposed of by drain.

- STORAGE -

Store concentrate in sealed container at ambient temperature. Do not freeze.

- PRECAUTIONARY STATEMENTS -

KEEP OUT OF REACH OF CHILDREN CAUTION

Contrex CA contains organic acid. Contents may cause eye irritation. Avoid contact with eyes.

-FIRST AID-

If in eyes: flush eyes with copious amounts of water for 15 minutes. Seek medical attention if eye irritation persists.

- PACKAGING -

1 Gallon /4 per case

- CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
1 Gallon	5701

For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com

®Decon Labs, Inc.

Revised 4/02/2012