



# Surface Disinfectant Wipes

## Technical Data Sheet

### EPA Reg. No. 8383-7-56753

**- BRIEF DESCRIPTION -**

*CiDecon Plus Wipes are disinfectant wipes in a convenient pop-up canister. An effective disinfectant and active deodorant for laboratories, hospitals, clinics, medical, dental and veterinary offices, industrial clean rooms, homes, nursing homes, ambulances, etc. For use on hard, non-porous, inanimate surfaces, they provide broad range efficacy against bacteria, virus and fungi. Proven effective against tuberculosis, HIV-1 and many other pathogenic organisms. CiDecon Plus Wipes are safe, effective and meet OSHA bloodborne pathogen standard.*

**- DIRECTIONS FOR DISPENSING -**

Pull up tab and remove cover from container. and pull up the wipe corner from the center. Twist the corner into a point and insert about one inch of the wipe through the X hole on the underside of the cover. Pull the wipe completely through and replace cover onto container. Pull out first wipe at a 90° angle. Remaining wipes feed automatically. Always reseal the hinged cap after each use to ensure moisture retention.

**- USE PROTOCOL -**

**To Disinfect and deodorize:** PRECLEAN surface to remove soil and filth. Thoroughly wet pre-cleaned surface and allow surface to remain wet for 10 minutes at room temperature to kill listed organisms.

**- ACTIVE INGREDIENTS -**

Phenol.....1.56 %  
 Sodium Phenate.....0.06 %  
 Inert Ingredients.....98.38 %

**- EFFICACY -**

PASSES AOAC EFFICACY STANDARDS FOR INSTITUTIONAL AND HOSPITAL DISINFECTION

- **Virucidal :** *Avian Influenza A virus (H9N2), , Influenza A<sub>2</sub> virus (Japan 305/57 Asian strain), Cytomegalovirus, Coronavirus, Polio type 1 virus, Vaccinia virus, Herpes simplex types 1/F and 2/G (oral, ocular and genital), HIV-1 (AIDS virus.)*
- **Bactericidal:** *Staphylococcus aureus, Salmonella enterica, , Pseudomonas aeruginosa, Methicillin resistant Staphylococcus aureus, Vancomycin resistant Enterococcus faecium (VRE), E. coli, Streptococcus pyogenes, Streptococcus salivarius, and Mycobacterium tuberculosis.*
- **Fungicidal:** *Trichophyton mentagrophytes (Athlete's Foot fungus).*

**- STORAGE AND DISPOSAL -**

Store at room temperature. Non- refillable container. Do not use or refill this container. Discard empty container in trash or offer for recycling if available.

# CiDecon® Plus Wipes Surface Disinfectant Wipes Technical Data Sheet

**FOR HOSPITAL DISINFECTION:** This product is not to be used as a terminal sterilant/high level disinfectant on any surface or instrument that (1) is introduced directly into the human body, either into or in contact with the bloodstream or normally sterile areas of the body, or (2) contacts intact mucous membranes but which does not ordinarily penetrate the blood barrier or otherwise enter normally sterile areas of the body. This product may be used to preclean or decontaminate critical or semi-critical medical devices prior to sterilization or high level disinfection.

**- PACKAGING -**

**180 wipes per canister / 12 canisters per case**

**- CATALOG NUMBER -**

| <u>Manufactured for:</u> | <u>Item</u> | <u>Cat. No.</u> |
|--------------------------|-------------|-----------------|
| Decon Labs, Inc.         | 180 wipes   | 8511            |

**- PRECAUTIONARY STATEMENT -**

**HAZARDS TO HUMANS AND DOMESTIC ANIMALS. KEEP OUT OF REACH OF CHILDREN. CAUTION:** Causes moderate eye irritation. Avoid contact with eyes or clothing. Wash thoroughly with soap and water after handling.

**FIRST AID:** If in eyes, remove contact lens. Hold eye open and slowly rinse with water. If irritation persists for more than 5 minutes continue rinsing and consult a physician or poison control center.

---

**For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

---

®Decon Labs, Inc. Revised 8/18/2011

**Page 2 of 2**