

DNase Displace™

Technical Data Sheet

- BRIEF DESCRIPTION -

DNase Displace removes DNase, and unwanted DNA. It works instantly so there is no waiting. DNase Displace eliminates the time-consuming method of autoclaving. DNase Displace can be used on all surfaces including glass, plastic, metal, and ceramic. DNase Displace is phosphate-free, safely disposable, non-toxic, and totally rinsable for the decontaminated surface.

- USE PROCEDURE -

Use DNase Displace full strength.

Method 1: Thoroughly wet item to be decontaminated with DNase Displace - ensure all surfaces are coated. Swirl for a few seconds, drain, and immediately rinse twice with the DNase-free water.

Method 2: Soak item in DNase Displace until needed. Drain and rinse as above.

Method 3: Soak an DNase-free wipe in DNase Displace, wipe over surface to be decontaminated. Rinse as above.

For total radioactive decontamination, soak item in DNase Displace for 2-12 hours. Rinse as above. Wipe over benchtops and sinks to remove residual contamination.

- CAUTION -

Alkaline! Severe eye irritant, mild skin irritant. Wear eye and skin protection when handling DNase Displace. In case of contact with eyes, immediately flush with water for fifteen minutes and contact a physician immediately.

- PACKAGING -

250 mL (8 oz.) / 12 per case

- CATALOG NUMBERS -

<u>Item</u>	<u>Fisher Cat. No.</u>	<u>FoxWest Sales Cat. No.</u>
8 oz.	04-355-135	1409

**For Additional Technical Information call
FoxWest Sales Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

FoxWest Sales, Inc.

Revised 12/05/2014