



Dri-Clean

Powdered Detergent



Dri-Clean™ is a powdered detergent for use in cleaning all labware and equipment. Use **Dri-Clean™** for soak, manual, and ultrasonic cleaning applications. Removes protein, dried blood, and most organic and inorganic contaminants from glassware, plasticware, ceramics, and metal surfaces. Completely free-rinsing and residue-free. **Dri-Clean™** is packaged in a wide-mouth plastic container with a convenient handle. This makes it easier to handle and prevents the absorption of water and caking associated with powdered detergents packaged in milk cartons.

WORKING SOLUTIONS: Dissolve 1/2 oz. (1 Tbsp.) in one gallon of water. For tenacious contaminants, use 1 oz. (2 Tbsp.) or more and dissolve in hot water. Clean by manually scrubbing, soaking, or use ultrasonic bath. For critical applications, clean and rinse with distilled or deionized water.

STORAGE: Store in original container with lid tightly sealed. Contamination with water or high humidity may cause caking.

PRECAUTIONARY STATEMENT:

Contains sodium carbonate. May cause eye and skin irritation. Avoid eye contact or prolonged skin contact. Avoid breathing dust.

WEAR PROTECTIVE EYEWARE AND GLOVES. KEEP OUT OF REACH OF CHILDREN

FIRST AID:

Eye Contact: Flush eyes with water for 15 minutes. Get medical attention if irritation persists.

Skin Contact: Rinse with water, remove contaminated clothing.

Ingestion: Give water or milk. Never give anything by mouth to an unconscious person. Get medical attention immediately.

Packaging

Size	Case QTY	Decon #
2kG (4.4lbs.)	6	2502
50kG (110lbs.)	-	2501

Rev. 12/2022