



PROGUARD®

Professional Handcream Technical Data Sheet

- BRIEF DESCRIPTION -

Proguard Professional Handcream combines emollients to moisturize dry, cracked skin with a barrier film that seals in the moisture, seals out the problems. It is particularly effective in eliminating the irritation caused by wearing latex gloves. Because Proguard does not contain perfumes, oils, greases, petroleum, aloe or cocoa butter, it is completely compatible with and will not compromise the protection of latex gloves.

- USE PROCEDURE -

Formulated Specifically For The Hands of Laboratory and Healthcare Professionals.

Most adverse/dry skin conditions will respond well to regular application of Proguard, and we recommend its use immediately after hand washing and always before gloving. Chronic skin disorders should be attended to by a physician. In the unlikely event of any skin reaction to Proguard, its use should be discontinued immediately.

Wash hands and dry. Apply one shot approximately 1 gram Proguard from the Dispenser Pump or a sparing amount from the tube to the palm of one hand. Rub in with a washing action to cover entire skin surface. Extremely dry skin may require a second application. Proguard absorbs rapidly to leave a silky, dry film. For best results, it should be used after every hand scrubbing and always before wearing protective gloves.

- PROGUARD COMPATIBILITY WITH LATEX GLOVES -

“Information presented by Dr. T. Lowe of the Food and Drug Administration (FDA) at the Fourth Conference on Occupational Hazards to Health Care Workers (University of Washington, May 1998) addressed not only the FDA’s current efforts to establish efficacy criteria for latex surgical and examination gloves, but also noted the significant deterioration of such gloves when exposed to petroleum-based lubricants...”

“...Therefore when hand creams are utilized, particular attention should be directed to the cream’s formulation and the effect it has on the barrier properties of the gloves.”

Federal Register, Volume 56;235, 12/6/91

Proguard Professional Handcream contains no petroleum products and no ingredients that deteriorate latex. If you wish to review our Latex Compatibility Test Procedures, call (800) 332-6647 and request a copy.

- PACKAGING -

3 oz. (85 g) Flip-Top Tube / 12 per case
16 oz. (450 g) Pump-Pack / 6 per case
Wall Mount Bracket for 16 oz. / 12 per case

- CATALOG NUMBERS -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	3 oz. (85 g)	9503
	16 oz. (450 g)	9516
	Wall Bracket	9517

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

®Decon Labs, Inc.

Revised 10/12/2011