



The Pill™

Autoclave Deodorizer Capsule

Technical Data Sheet

- BRIEF DESCRIPTION -

THE PILL Autoclave Deodorizer Capsule neutralizes odors; it is not simply a perfumed mask. Each PILL contains concentrated odor neutralizer liquid in a gelatin capsule that keeps pungent deodorizer fluids off hands and clothes and makes for accurate and economical unit dosing. One PILL deodorizes seven cubic feet of autoclave capacity with full load of soiled labware. (Not suited to small tabletop autoclaves).

- USER PROTOCOL -

Use one PILL per 7 cubic ft. autoclave capacity. Place in a beaker with one-inch cold water. Place beaker in bottom of autoclave with soiled load and run cycle.

For non-autoclave use, e.g. deodorizing animal rooms, drop one PILL into large beaker or bucket full of very hot water. Allow to vaporize. If possible, circulate air by way of a fan.

- CONTRA-INDICATIONS -

One PILL is too strong for small tabletop autoclaves. There will be a pungent residual odor of unused deodorizer if used in units much smaller than 5 cubic ft. capacity.

- PACKAGING -

100 capsules / 20 containers per case
500 capsules / 6 containers per case

- CATALOG NUMBERS -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	100 Capsules	5100
	500 Capsules	5500

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

™ Decon Labs, Inc. Revised 11/2006