



CITRAD™

Acid Cleaner & Detergent



Citrad™ concentrated acidic detergent is ideal for the removal of salts, scale, and metal-oxide. Also used for metal brightening and pharmaceutical-process equipment. Citrad™ contains a safe, organic acid and is phosphate-free, so it will not interfere with sensitive equipment or testing procedures. Citrad™ is also completely free-rinsing and biodegradable.

Citrad™ concentrate must be diluted with distilled, deionized or clean potable water and can be used as a soak, scrub or ultrasonic cleaner.

Citrad™ is used to clean labware, stills, cleanrooms, optical parts, pharmaceutical equipment, and manufacturing equipment. Citrad™ is also effective in removing metal oxides, scale, soil, grease, chemicals and solvents and other compounds from surfaces and equipment.

DIRECTIONS FOR USE: Add 1.25 ounces to 2.5 ounces per gallon of water (1% to 2% v/v). For more difficult applications increase the concentration of Citrad™ solution and/or increase the temperature.

Citrad™ will foam if dispensed by sprayer.

DISPOSAL: All ingredients in Citrad™ are biodegradable and are considered safe when recommended use dilutions are disposed of by drain.

**1 Gallon
4 per case
Decon Catalog No. 4401**

Rev. 11/2022