



Contrex® AP is a powdered labware detergent. It removes virtually any contaminant including grease, oils, blood, tissue, particulates, deposits, chemicals and solvents. **Contrex® AP** works by soaking, scrubbing, or in ultrasonic cleaning baths. **Contrex AP** is totally rinsable eliminating detergent build-up that interferes with analysis, tissue culture, and other critical procedures. Its activity can be accelerated by heating, agitation, and ultrasonics, or by increasing concentrations. **Contrex® AP** is biodegradable and readily disposable.

Packaging		
Size	Case QTY	Decon #
4 lbs.	9	5204
25 lbs.	-	5225

APPLICATIONS: **Contrex® AP** is not to be used with machines, product will foam. After cleaning, rinse thoroughly. Safe on most surfaces including glass, plastic, and most metals. Cleaning applications include:

- **Labware**
- **Instruments**
- **Equipment**
- **Production Mixers, fillers, storage containers**
- **Optics**

WORKING SOLUTIONS: Make a fresh 1% **Contrex® AP** working solution (1.25oz. per gallon) in deionized or distilled water for best results. Soak, scrub or sonicate as needed.

RINSABILITY: **Contrex® AP** is totally rinsable - however, numerous factors affect rinsability. These factors include water quality, temperature, volume of water, nature of surface, shape/size of the vessel.

STORAGE: Keep contents in a clean and dry environment.



For details on Decon's Full Line of Contrex Detergents visit



Rev. 09/2022