



Contrex® AW is a low-foaming, concentrated, powdered autowasher detergent for use in all automatic labware washing machines. It is extremely effective at removing virtually any contaminant from the surface of labware, scientific apparatus, and process equipment. **Contrex® AW** is free-rinsing - eliminating detergent build-up in the automatic machines.

APPLICATIONS: When using **Contrex® AW** always follow machine manufacturer's instructions. Typical cleaning applications include:

- **Labware**
- **Instruments**
- **Process equipment**
- **Glass & ceramics**
- **Optics**

WORKING SOLUTIONS: Typical use is approximately 1.25 ounces of solution per gallon of water (1% solutions) used during the wash cycle. For more difficult applications increase the concentration and/or temperature. Heavily soiled labware may need a prewash using **Contrex® AL**. Ensure adequate rinse cycle for spot-free finish. All cleaners work best in deionized or distilled water.

RINSABILITY: **Contrex® AW** is totally rinsable - however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water.

STORAGE: Keep contents in a clean and dry environment.

Packaging

Size	Case QTY	Decon #
4 lbs.	9	5604
25 lbs.	-	5625
50 lbs.	-	5650



For details on Decon's Full Line of Contrex Detergents visit

www.deconlabs.com

Rev. 09/2022