



Contrex® CA concentrated acidic detergent is formulated for use in descaling, mineral removal, oxide removal, or in applications where an acidic detergent is indicated. Contrex® CA contains a safe, organic acid and is phosphate-free, so it will not interfere with sensitive equipment or testing procedures. Contrex® CA is also completely free-rinsing and biodegradable.

**Decon Catalog Number
5701
4 x 1 Gallon**

APPLICATIONS: Contrex® CA concentrate can be diluted with distilled, deionized, or clean potable water and can be used as a soak, scrub, or ultrasonic cleaner.

Contrex® CA is not for use in machines - product will foam. In order to accelerate cleaning - increase the solution strength or temperature. Contrex® CA is safe on most surfaces including glass, plastic, and most metals. Cleaning applications include:

- **Labware**
- **Instruments**
- **Equipment**
- **Production Mixers, fillers, storage containers**
- **Optics**

WORKING SOLUTIONS: Make a fresh 1 to 2% Contrex® CA working solution (1.25 to 2.5 oz. per gallon) in deionized or distilled water for best results. Clean by scrubbing, soaking, or by ultrasonic cleaning. For critical applications use deionized water. After cleaning, rinse thoroughly.

RINSABILITY: Contrex® CA is totally rinsable - however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape and size of the vessel.

DISPOSAL: All ingredients in Contrex® CA are biodegradable and are considered safe when recommended use-dilutions are disposed of by drain.

STORAGE: Store concentrate in sealed container at ambient temperature. Do not freeze.



**For details on Decon's Full Line
of Contrex Detergents visit**

www.deconlabs.com

Rev. 09/2022