



Technical Data Sheet

- BRIEF DESCRIPTION -

Contrex CF is a cation free and a phosphate free liquid detergent concentrate. Contrex CF is formulated for use in cleaning of equipment and electronics that are sensitive to sodium and metal cations and also phosphates.

- APPLICATIONS -

Contrex CF is totally rinsable eliminating detergent build-up that interferes with analysis and other critical procedures.

To accelerate cleaning increase concentration and/or temperature. After cleaning, rinse thoroughly. Cleaning applications include:

- Electronic components
- Solder/Flux removal
- Environmental lab ware

- WORKING SOLUTIONS -

Contrex CF may be diluted with distilled, deionized or clean potable water and can be used as a soak, scrub, ultrasonic cleaner, or machine washing. Make up a fresh 3 % to 5 % Contrex CF working solution (4 to 6 oz. per gallon) in deionized or distilled water for best results. Solutions of Contrex CF are biodegradable and readily disposable. Protective gloves and safety glasses are recommended.

- RINSABILITY -

Contrex CF is totally rinsable – however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape and size of the vessel.

-STORAGE -

Store concentrate in sealed container at ambient temperature. Do not freeze.

- PRECAUTIONARY STATEMENTS -

KEEP OUT OF REACH OF CHILDRREN

CAUTION

Contrex CF contains solvents. Can cause eye irritation. Avoid contact with eyes. Wear gloves, eye protection and avoid breathing vapors.

-FIRST AID-

If in eyes: flush eyes with copious amounts of water for 15 minutes. Seek medical attention if eye irritation persists.

- PACKAGING -

1 Gallon / 4 per case

- CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
1 Gallon	5801

For Additional Technical Information call
Decon Customer Service
 (800) 332-6647, or
 See our website at www.deconlabs.com

©Decon Labs, Inc.

Revised 4/02/2012