



Contrex® EZ is a concentrated, powdered enzymatic detergent. It is formulated for use as an enzymatic detergent for use in manual or soak cleaning of labware and equipment contaminated with protein and fats.

APPLICATIONS:

Contrex® EZ is not for use in machines
- product will foam.

Safe on most surfaces including glass, plastic, and most metals. Cleaning applications include:

- **Labware**
- **Instruments**
- **Endoscopes**
- **Medical Devices**
- **Process equipment**
- **Food & Dairy surfaces**
- **Ultra filtration & reverse osmosis filters**

Packaging

Size	Case QTY	Decon #
4 lbs.	9	5404
25 lbs.	-	5425

WORKING SOLUTIONS: Contrex® EZ concentrate can be diluted with distilled, deionized, or clean potable water. Make up a fresh 1% Contrex® EZ working solution (1.25oz. per Gallon) in deionized or distilled water for best results. Clean by scrubbing, soaking, or by ultrasonic cleaning.

RINSABILITY: Contrex® EZ is totally rinsable - however - numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape, and size of the vessel.

STORAGE: Keep contents in a clean and dry environment.



**For details on Decon's Full Line
of Contrex Detergents visit**

www.deconlabs.com

Rev. 09/2022