



OdorGone®

Autoclave Deodorizer Capsule

Technical Data Sheet

- BRIEF DESCRIPTION -

ODORGONE® Autoclave Deodorizer Capsules contain concentrated liquid odor neutralizers which absorb odors instead of just masking them. Gelatin capsule keeps deodorizing liquid off hands and clothes and makes for accurate unit dosing. Available in three pleasant fragrances.

- PACKAGING -

25 capsule pack / 8 packs per case
100 capsule pack / 6 packs per case

- USE PROTOCOL -

Place one ODORGONE capsule in a beaker and cover with one inch of cold water. Place beaker in bottom of autoclave with soiled load and run cycle. Alternately, you may place one ODORGONE capsule in each autoclave bag containing autoclaveable waste. Place bags in autoclave and run cycle.

- STORAGE -

Keep ODORGONE capsules sealed in original container. Store in a cool, dry place.

- CATALOG NUMBERS -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.		
	25 Capsules	
	Apple	APP25
	Lemon	LEM25
	100 Capsules	
	Apple	APP100
	Lemon	LEM100

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

®Decon Labs, Inc.

Revised 2/2008

-PRECAUTIONARY STATEMENT -

DO NOT INGEST. For use in autoclaves only.