



OdorGone®

Autoclave Deodorizer Capsule



ODORGONE® Autoclave Deodorizer Capsules contain concentrated liquid odor neutralizers which absorb odors instead of just masking them. Gelatin capsule keeps deodorizing liquid off hands and clothes and makes for accurate unit dosing. One capsule deodorizes 7 cubic feet of autoclave capacity with full load of soiled labware.

Note: Not suited to small tabletop autoclaves. One capsule is too strong for small tabletop autoclaves. A pungent residual odor of unused deodorizer will be present in units smaller than 5 cubic feet.

Available in two pleasant fragrances.

USE PROTOCOL: Use one **ODORGONE®** capsule per 7 cubic feet autoclave capacity. Place capsule(s) in a beaker and cover with one inch of cold water. Place beaker in bottom of autoclave with soiled load and run cycle. Alternately, you may place one **ODORGONE®** capsule in each autoclave bag containing autoclaveable waste. Place bags in autoclave and run cycle. For non-autoclave use, e.g. deodorizing animal rooms, drop one capsule into a large beaker or bucket full of very hot water. Allow to vaporize. If possible, circulate air by way of fan.

STORAGE & DISPOSOL: Keep **ODORGONE®** capsules sealed in original container. Store in a cool, dry place.

PRECAUTIONARY STATEMENTS: DO NOT INGEST. For use in autoclaves only.

PACKAGING

Scent	PACK QTY	CASE QTY	Decon Catalog #
Apple	25 Capsules	8	APP25
	100 Capsules	6	APP100
Lemon	25 Capsules	8	LEM25
	100 Capsules	6	LEM100

Rev. 12/2022