



SANIHOL[®] 70

70% Denatured Ethanol Solution



Sanihol[®] 70 is a 70% (v/v) denatured ethanol solution that is filtered to 0.2 microns. **Sanihol[®] 70** can be used to clean surfaces in microbiology labs and tissue culture labs. It is also effective in cleaning surfaces and process equipment in pharmaceutical, biotechnology, and medical device, cosmetic and production areas. The alcohol solution evaporates away completely so no rinse is required.

Use Protocol: Use in well-ventilated area. Spray or wipe solution on surface to be cleaned until it is completely covered. Allow to air dry. When using trigger spray, point away from eyes and skin. Hold bottle six to eight inches from surface to be treated.

INGREDIENTS

63.4% Ethyl Alcohol

3.2% Methyl Alcohol

3.4% Isopropyl Alcohol

30% USP Purified Water % by volume

Precautionary Statements

Flammable - do not come in contact with flame or extreme heat. Avoid eye and skin contact and ingestion.

SHELF LIFE
2 YEARS

Size	Case Qty	Decon Cat #
16 oz.	12	8616
1 Gallon	12	8601
55 Gallon	-	8655

Rev. 07/2021